

Method Abstract #77

Conductivity


Scope and Application

This method determines the conductance of water samples. This method conforms to Standard Method 2510 B, ASTM D1125, ISO 7888 and EN 27888

Method Summary

A conductivity cell consisting of two plates is submerged into a solution and a voltage is passed across the two plates. Ions in the solution are attracted to the plate of opposite charge and move between the two plates depending on the resistance of the solution. Conductivity is then expressed as the reciprocal of the electrical resistance and is commonly reported in units of microsiemens (μS). Prior to sample analysis, a calibration is performed using standard potassium chloride solution. The conductivity reading is then read directly from the conductivity meter.

Sample Calibration Curve

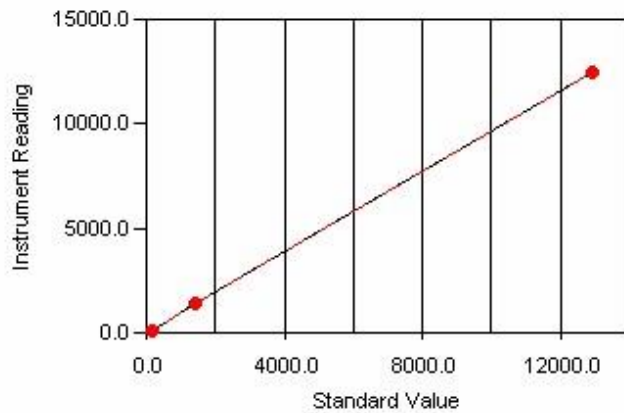

Method Performance

Parameter	Specification
Measuring Range*	0.01 – 199,999 μS
RSD for deionized water, 0.95 μS	19.10% or +/- 0.18 μS
RSD @ 147 μS	0.70% or +/- 1.03 μS
RSD @ 1413 μS	0.18% or +/- 2.54 μS
RSD @ 12,900 μS	0.11% or +/- 14.19 μS
RSD for Tap water, 939.90 μS	0.33% or +/- 3.10 μS

*Data for this measuring range was obtained using laboratory prepared standards formulated from potassium chloride. The measuring range may be increased by using auto-dilution

RSD Values are better than those specified in Standard Methods