



Instructions for Setting Up the MANTECH AutoMax 73 Autosampler

During shipment, a screw locks the horizontal arm into place.

1. Remove the black cap covering a screw hole on the right side of the sampler. Using a large Philips head screwdriver, remove the long screw, securing the y-arm to the right side of the sampler. Use the hole at the back of the autosampler to safely store the screw.

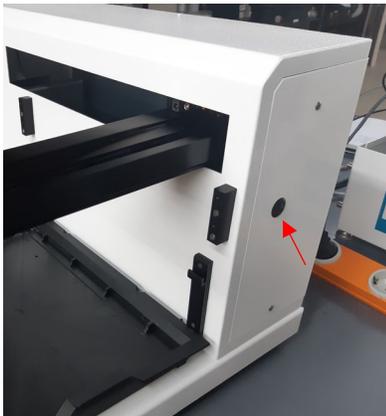

2. Place the black drip tray onto the base of the sampler.

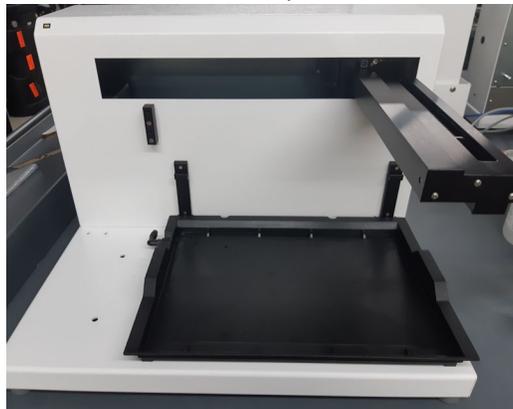


3. Plug the power and communication cables (PC-1205) into the back of the sampler.




4. Locate the z-arm assembly from the sampler box and two screws - if they are screwed into the bottom of the z-arm, remove them.



5. Remove the two right screws and loosen the far-left screw from the sampler arm.

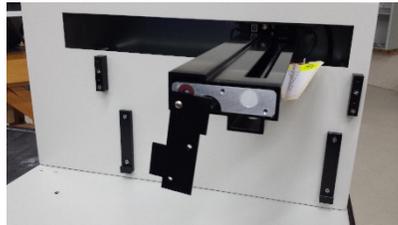

6. Slide the rod out to place the Z arm in and then slide the rod back in.

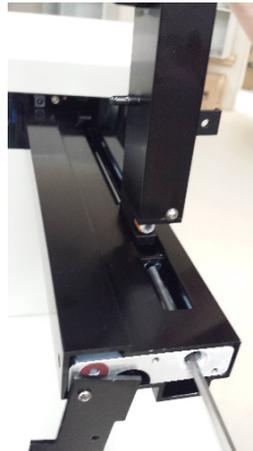


7. Screw the front plate back on.




8. Slide the Z arm towards the front of the arm to line up with the holes underneath and secure the Z arm with the two screws from Step 4.



9. Slide the z-arm back slightly to reveal the hole for the stability guide. Insert the stability guide into the slot and secure with the screw provided with the guide.

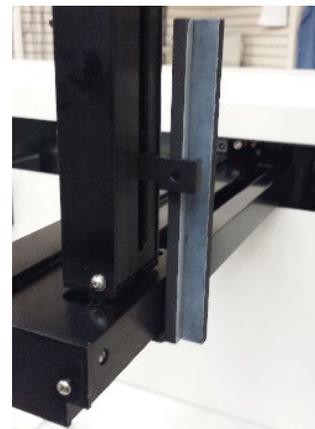

10. Attach the stiffener assembly.

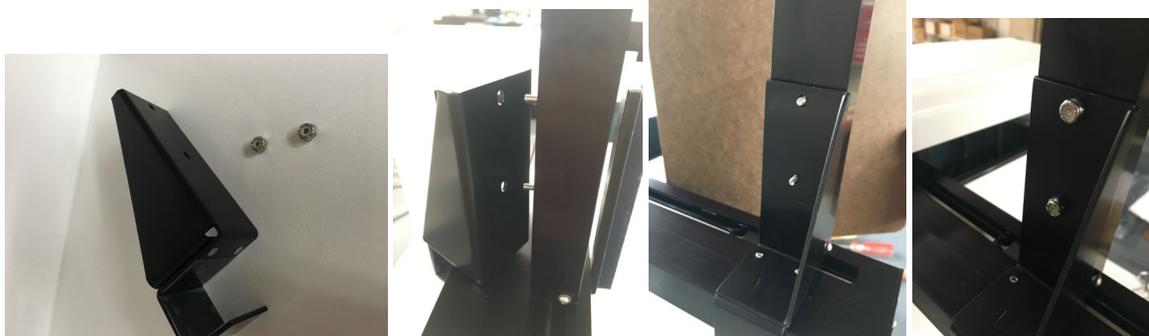

11. The AutoMax73 setup is now complete.

Note: Before packing the Autosampler for shipment, always secure the horizontal arm using the arm locking screw.



For any questions on setting up the AutoMax73 or other MANTECH materials, please contact your local distributor or: support@mantech-inc.com.

Document Change Log

Version	Date	Author	Changes
2	21-Aug-2019	Heather Jasumani	<ul style="list-style-type: none">Added step for stiffener assembly



www.mantech-inc.com

MANTECH Inc.
5473 Highway 6 North
Guelph, Ontario N1H 6J2
Canada
519-763-4245