

Shortcuts | Alert Configuration | Run Configuration

Calibration



New Sensor (Break-In)



QC Routine



Secondary Standards



Sample Analysis



Performs 3 Calibrations and a QC-Check (measures a COD value).

Performs 4 Calibrations and a QC-Check.
Used ideally for new sensors. (Generally, only needed to be used once).

Performs an Analysis and re-calibrates and analyze if the initial analysis failed.
Used for verifying if calibration is still valid.

Performs an analysis to check if secondary standard is within range of calibration.
Useful for verifying calibrations/QC Routines.

Performs any number of analyses based on what the user desires.

Notes:

- Number of shortcuts present on your database will be based on the COD range(s) purchased.
- For further questions/inquiries, please contact: support@mantech-inc.com