

IntelliPeri™ Estimated Flow Rates by Speed and Tubing Size

This document outlines the estimated flow rate of the MANTECH IntelliPeri™ for various speed settings and tubing diameters. For more information please contact us at support@mantech-inc.com.

Tubing Diameter (inches)	Speed Setting (%)	Flow Rate (mL/min)
1/8"	6 (min)	1.9
	10	17.88
	20	56.22
	30	92.27
	40	126.02
	50	157.47
	60	186.63
	70	213.49
	80	239.05
	90	260.32
	98 (max)	276.48
1/4"	3 (min)	2.79
	10	61.82
	20	146.15
	30	230.48
	40	314.81
	50	399.14
	60	483.47
	70	567.80
	80	652.13
	90	736.46
	99 (max)	812.36