



# MANTECH

OPTIMIZE YOUR RESULTS. PROTECT OUR ENVIRONMENT.




## AM500 SERIES ANALYZER

FOR METALS DIGESTION SCREENING & SAMPLE SPLITTING

# INCREASE AUTOMATION, REDUCE AILMENTS FROM MANUAL TASKS

MANTECH's AM500 series analyzers take automation to the next level by eliminating tedious, repetitive aliquoting and analysis steps required for metals digestion screening & other routine analyses.

WATCH SYSTEM IN ACTION

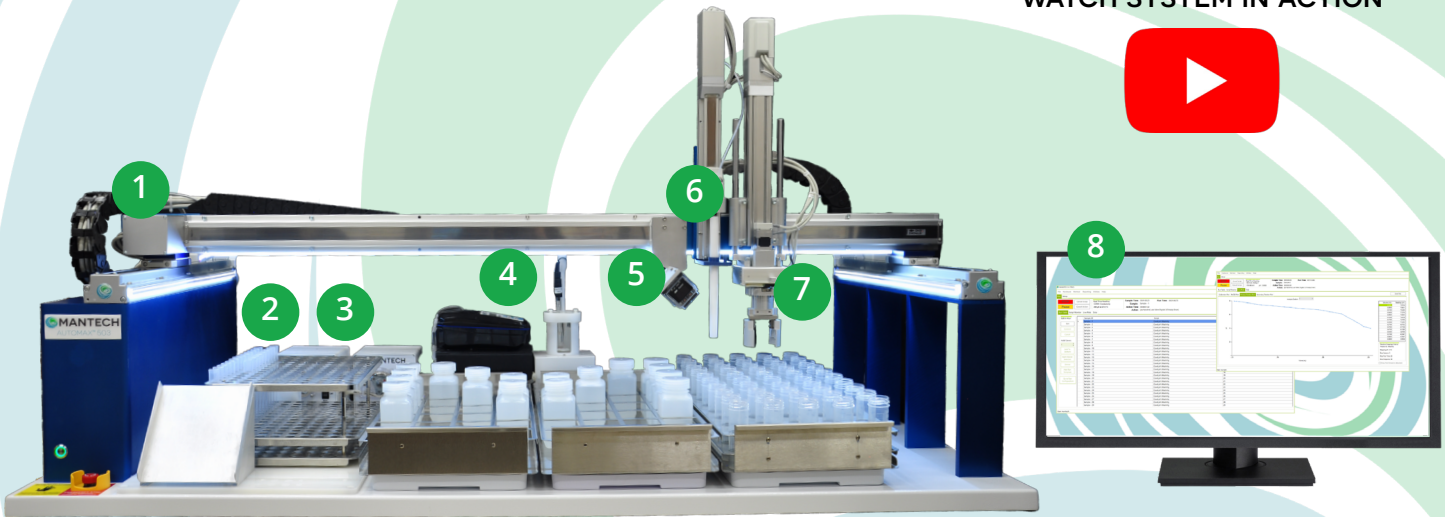

## AM500 SERIES ANALYZER FEATURES

- 1 H-Frame Autosampler**  
Includes two z-arms for analysis. Maintenance free with fully sealed, self-lubricating axes.
- 2 IntelliRinse Station**  
Cleans all wetted analysis vessels and sample lines with fresh water before proceeding to next sample to eliminate cross-contamination.
- 3 Intelligent Pumps**  
Used to deliver fresh water for rinsing and aliquots sample into external vessels. Communicate directly with software.
- 4 External Analysis Vessels**  
Analysis occurs within a TitraSip™ vessel and a T10 turbidity meter so the remaining sample can be used for other tests.
- 5 Barcode Scanner**  
Scans barcodes regardless of position on bottle (upside down, sideways, etc.). Pulls key information from LIMS. Ensures correct pipette tip, bottle or lid is picked up.
- 6 Pipette Tip Pickup & Ejector**  
Disposable or reusable pipettes available. Built-in sensor ensures pipette has been picked up or ejected. Pipette is ejected into an external disposal bin after analysis.
- 7 Rotating Gripper**  
Grips and rotates lids to cap and uncap field bottles. Built-in sensor ensures cap is removed or placed on.
- 8 MANTech Pro Software**  
User-friendly interface for reviewing current and historical data, generating reports, sending email alerts and more.

**AUTOMATE ADDITIONAL PARAMETERS  
CONDUCTIVITY, FLUORIDE, ORP, SULFIDE AND MORE!**

Request a quote: [info@mantech-inc.com](mailto:info@mantech-inc.com)

# Metals Digestion Screening Cycle Time ~ 3 Minutes



In this time, the AM500 series analyzer:


- Scans barcodes
- Unscrews caps
- Transfers sample to external analysis vessels
- Measures pH and turbidity
- Adds HNO<sub>3</sub> if needed
- Performs cleaning procedures
- Records results in software

**Add Sample Splitting and the Total Cycle Time ~ 7 minutes**

## ANALYSIS COMPONENTS



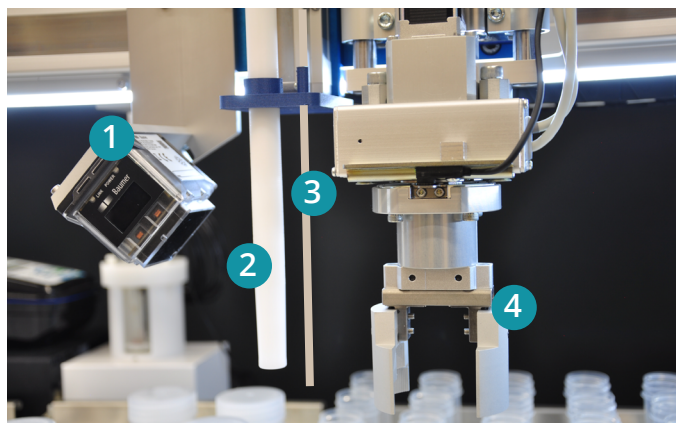


The AM500 series analyzers includes intelligent modules that communicate bi-directionally with MANTECH Pro software.

### ARM COMPONENTS



1. **Barcode Scanner** scans barcodes and ensures pipette tips, bottles and lids are picked up correctly.
2. **Pipette Pickup and Ejector** picks up a unique pipette tip and ejects it after analysis.
3. **Reagent Dispense Tip** dispenses HNO<sub>3</sub> into sample as needed.
4. **Rotating Gripper** uncaps and caps field bottle lids.


### PUMPS & EXTERNAL ANALYSIS VESSELS

				
<b>IntelliPeri™ Pumps</b> delivers rinse water and drains vessels.	<b>IntelliDose™ Pump</b> Class A pipetting for aliquoting samples.	<b>100,000 Step Buret</b> delivers precise volume of reagents.	<b>TitraSip™ Vessel</b> houses pH electrode & stirrer for analysis.	<b>T10 Turbidity Meter</b> measures turbidity with 30mL of sample.

Request a quote: [info@mantech-inc.com](mailto:info@mantech-inc.com)

# SOFTWARE & EXTERNAL COMPONENTS

## MANTECH PRO SOFTWARE

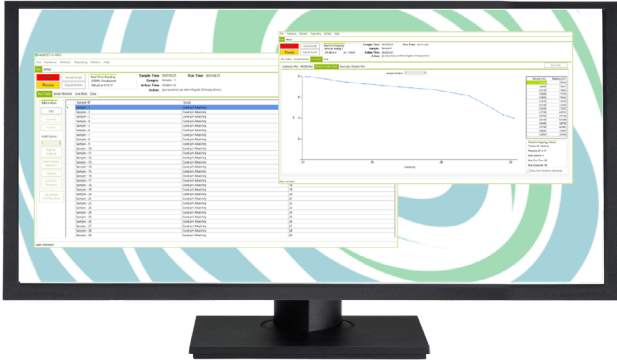

- Built for Windows 10 & 11 Pro OS
- Allows auto-security updates
- Shortcut buttons
- Import/export capability to LIMS
- Live view of data
- Historical data search
- Email alerts
- User account permissions
- Context sensitive Help Menu


## EXTERNAL COMPONENT OPTIONS

**Safety Shield with Safety Interlock** shields the exterior of the autosampler.

**Signal Light** includes three colors. The signification of each color can be customized to indicate system ready for analysis, run in progress, run ended, etc.

**LED Light Strips** illuminate the autosampler.


**Table with Casters** includes storage space on lower shelf.

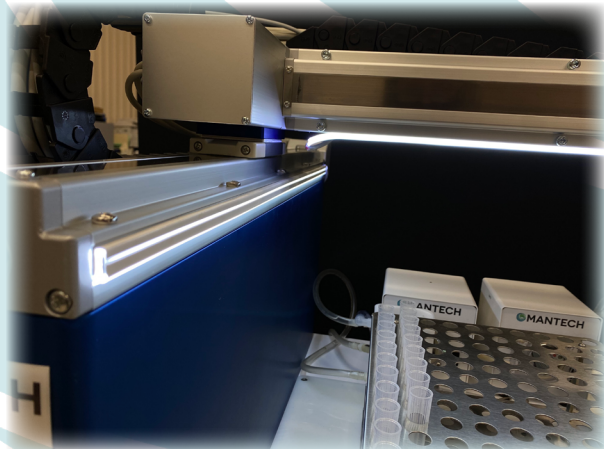



Request a quote: [info@mantech-inc.com](mailto:info@mantech-inc.com)

## AM500 SERIES DIMENSIONS

There are two autosampler sizes available for the AM500 series, differing in sample rack capacity, the AM503 can hold 3 sample racks while the AM505 can hold 5.

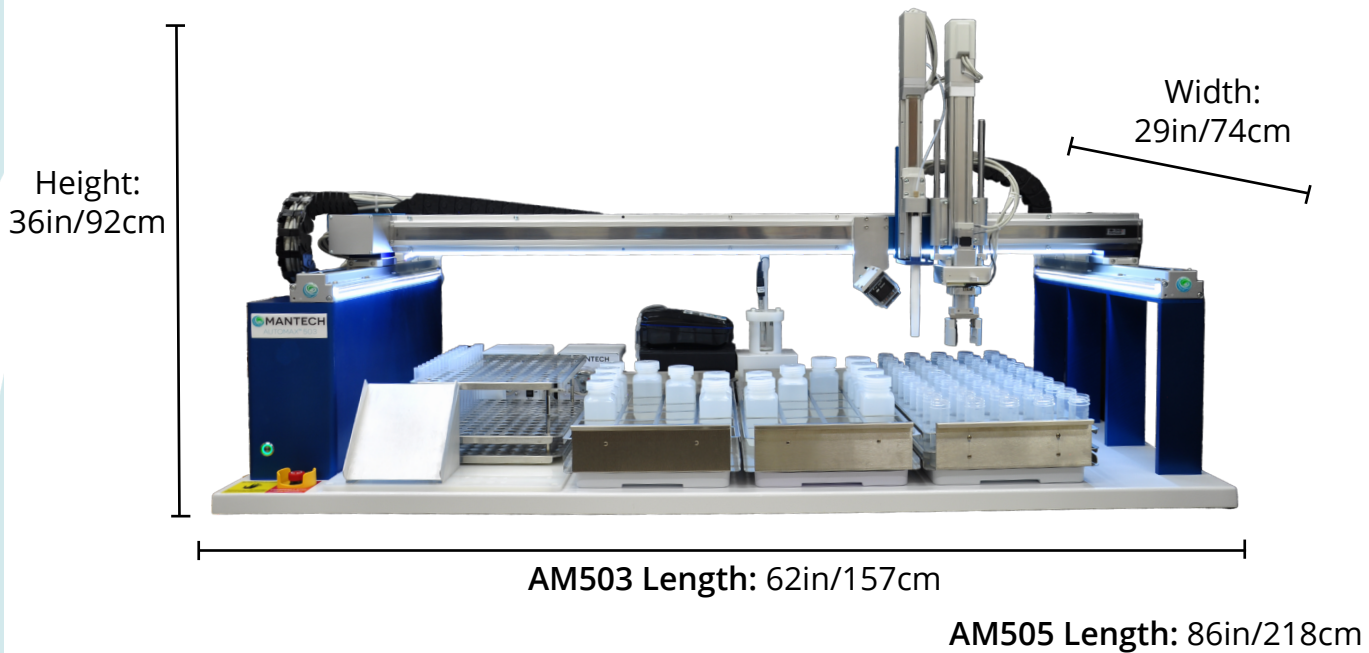

*Note: Enclosure adds 4" to the length and width of the autosampler.*

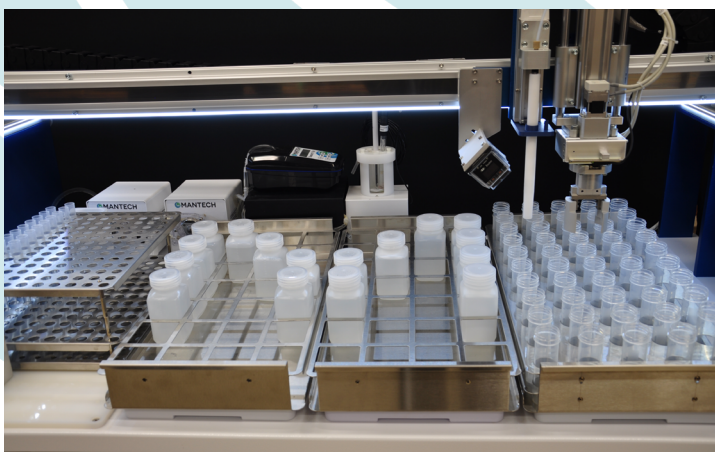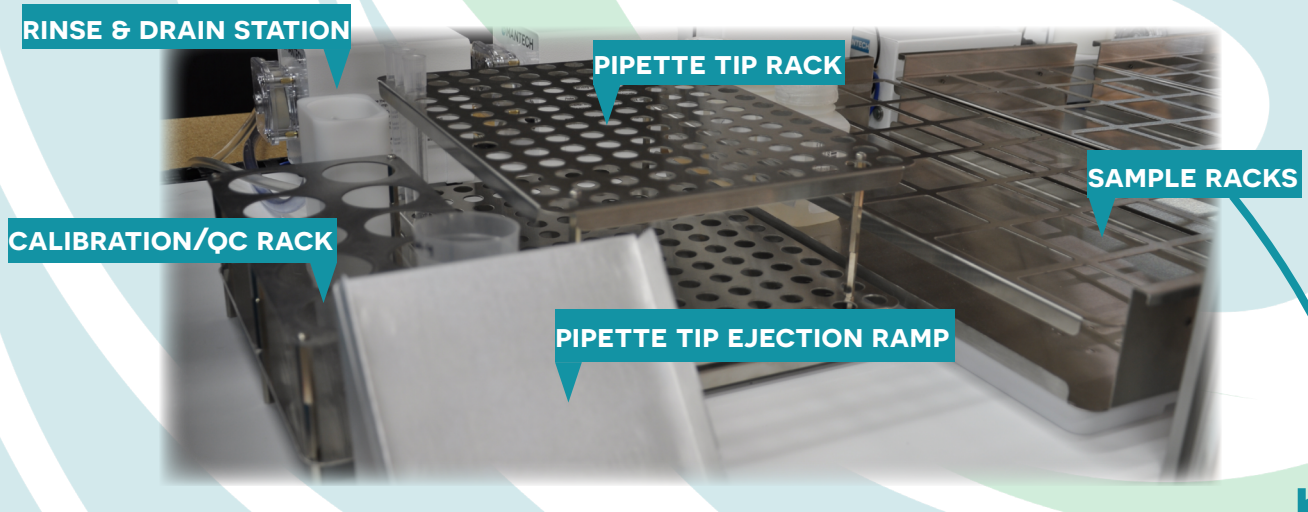

CAPACITIES PER RACK	
1L Bottles	14 positions
500mL Bottles	20 positions
250mL Bottles	28 positions
50mL Tubes	55 positions
15mL Tubes	242 positions

Request a quote: [info@mantech-inc.com](mailto:info@mantech-inc.com)